
Operating Manual

Control SM 7



Luigs & Neumann GmbH
Boschstrasse 19 * D-40880 Ratingen * Germany
Tel. : +49 2102 947000 * Fax.: +49 2102 442036
info@luigs-neumann.com * www.luigs-neumann.com

Control SM7

The control SM7 is operated with the keypad SM7 and is very versatile. Depending on the assembly, 1 to 9 motor-driven LN-axes can be connected.

Three slots are available for inserting circuit boards in a control box.

There are three different circuit boards:

Circuit boards with one, two or three motor-drivers.

Free slots can be retrofitted with circuit boards.

The control SM7 is also compatible with the control systems SM5, SM6 and SM8.

Up to eight LN-controls can be networked in-line.

The control SM7 features 3 operating modes:

A: The control SM7 is used in combination with one keypad.

B: The control SM7 is used in combination with two keypads.

C: The control SM7 is integrated into an existing SM5/SM6/SM7/SM8 system.

Contents

Page

Introduction / Contents.....	1
Safety instructions.....	2
Overview: Control SM7.....	3
Start up of the control SM7	4 / 5
Implement control SM7 into an existing LN control system.....	6
LN-SM7 ConfigurationTool.....	7
Changing address.....	7
Setting the control address.....	8 / 9
Example.....	10
Example: Address the Slave-control on 114.....	11 / 12
Example: Configuration motor connections.....	13 / 14 / 15 / 16
Appendix A, RS232/V.24.....	17
Technical data.....	18

Safety instructions

1.1 General instructions

The manipulators and moving-tables are designated for the positioning of repositories, microscopes or tools. (e.g.: capillaries, measuring electrodes, stimulating electrodes... etc.)

For a safe function of the manipulators pay attention to the operating and assembly manuals. Our service team will assist you, if additional information is required.

Comply with the security advice of this manual.

The intended protection can be endangered if the device is not used according to the operating manual of the producer.

1.2 User instructions

Do not touch the motor-driven manipulators during the positioning process in order to avoid injuries and bruises, and to avoid damaging the stated functions of the device.

As a result of the modular, individually constructed device by the customer, different areas can be danger zones with an increased risk of injury or wounds.

Avoid bringing your face too close to the moving devices, because the cramped space in combination with faulty operation of the manipulators can easily lead to the breakage of glass and glass splinters in the eye. The provided protective covering must be assembled according to the operation manual.

1.3 Transport instructions

The manipulators and moving tables are transported in a special packing in order to avoid possible damages during transportation.

1.4 Service and maintenance instructions

The manipulators and moving tables are maintenance-free.

In order to maintain the functional efficiency of the the devices, they must be protected from humidity and excessive heat. Strong jerks can lead to an incorrect adjustment which limits the functional efficiency, with the exception of the activities mentioned in the operating manual or if instructed by our service team. No personal changes are allowed.

1.5 Installation location instructions

Install the devices at locations with adequate air supply for aeration of the equipment.

The main control switch of the device has to be easily accessible at all times.

1.6 Disposal instructions

Broken devices or those no longer needed do not belong in the household rubbish! Dispose of them according to the local legal regulations. If in doubt, ask the service team of Luigs & Neumann for help.


1.7 Guarantee instructions

The producers are not liable for damages caused by unauthorized interference. Unauthorized interference terminates all warranty claims.

1.8 Repair and readjustment instructions


The devices must be packed according to the legal regulations. Contaminated devices must be cleaned beforehand if possible, otherwise a user declaration must be provided in which the substance is described and the absolute safety of the health of human-beings is confirmed.

Overview: Control SM7

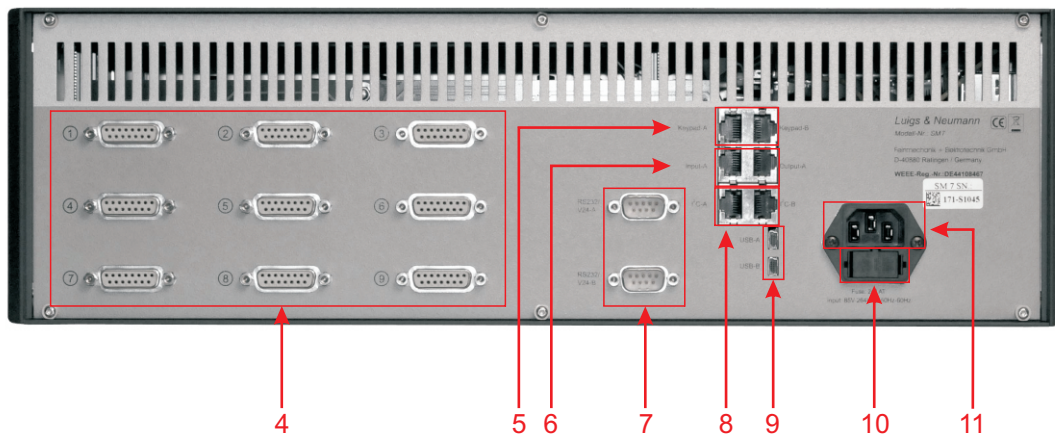


- 1 **Power switch**..... : On / off switch of the control
- 2 **Earth plug socket**..... : Earth the control box
- 3 **Display** : Display
- 4 **Motor connection 1 - 9** : Connection for motor-driven LN-axes
- 5 **Keypad A / B** : Plug for keypad SM7
- 6 **Input / Output** : Link SM7 with other controls
- 7 **RS232 / V.24** : Interface control
- 8 **I²C-BUS** : Add-on for additional devices
- 9 **USB** : Configuration of the controller and Interface control
- 10 **Fuse insertion**..... : Fuses: 2x T2A
- 11 **Power input**..... : 90-240VAC

Start up of the control SM7

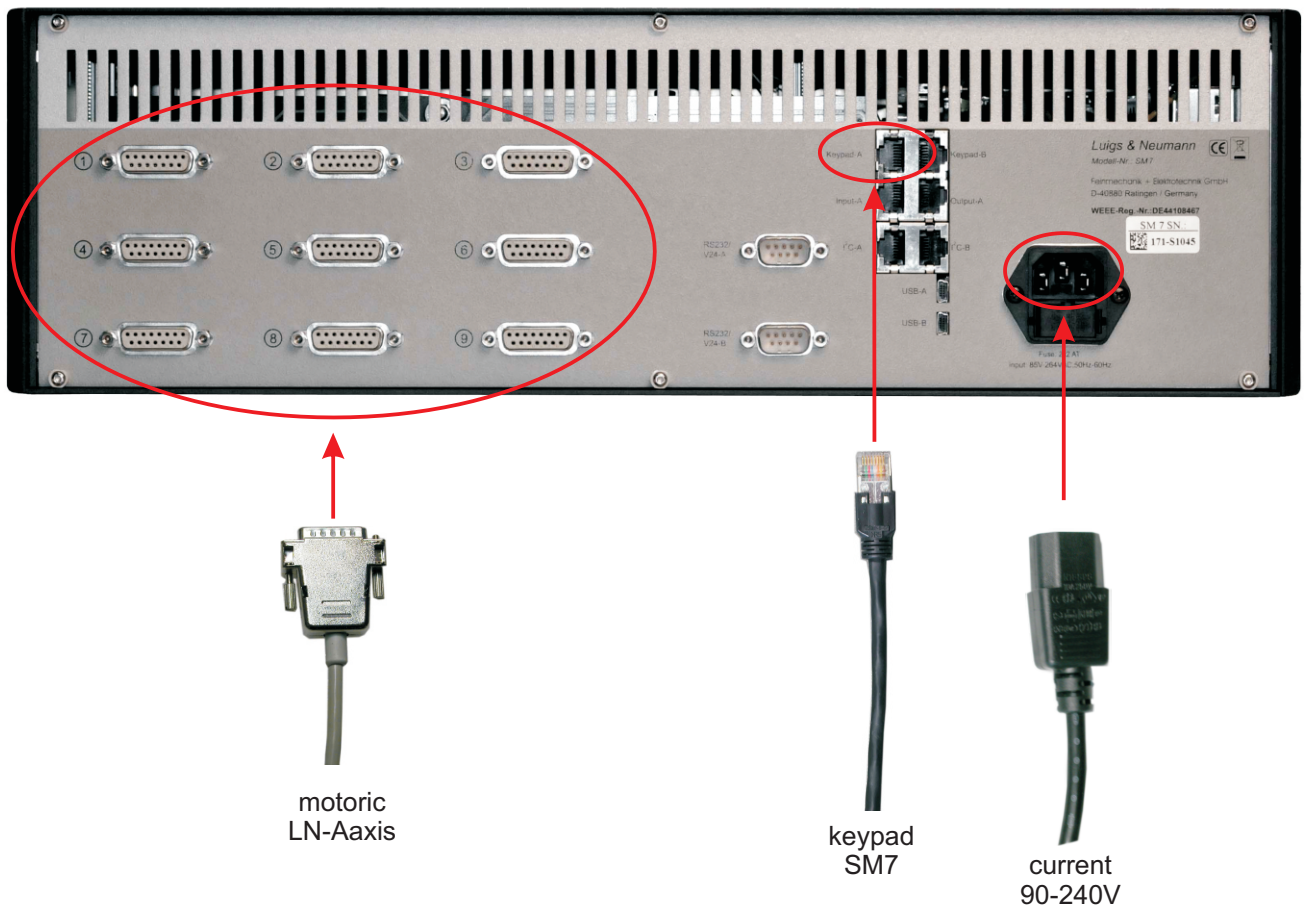

Connect the keypad and manipulators, only in the off-mode.
The keypad SM7 can only be connected to a control box with the adress 113.

- * Connect the control SM7 with the power cord.
- * Connect keypad SM7 with the keypad cable on the socket [5] Keypad-A. When using two keypads, connect the second keypad with the socket keypad B.
- * Connect the motoric LN-axes with the corresponding motor connections ①...⑨ of control SM7. The motor connections and the LN axes are colour-coded.

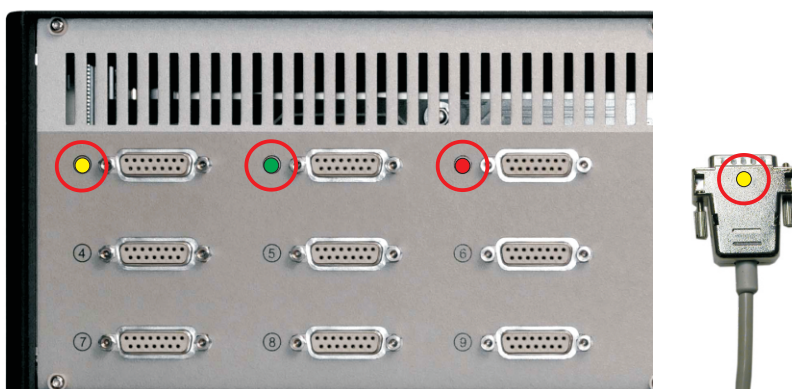

Start up of the control SM7

* Switch on the control SM7 over the power switch [1].

The display shows:

S	S	K		V	e	r	.	:	0	.	1	.	4			
S	e	a	r	c	h	i	n	g		s	t	a	g	e	s	

The program version of the Interface card.

S	y	s	t	e	m		r	e	a	d	y					
A	d	r	:	1	1	3	-	M	a	s	t	e	r			

The control address is shown.

The control is ready for operation.

Implement control SM7 into an existing LN control system

The control SM7 can be cross-linked with other controls (SM5, SM6, SM7 and SM8).

Up to 8 controls can be linked which can be operated with the keypad SM7.

The first control which is connected with the keypad must have the address 113 and is labelled as Master control.

All subsequent additional controllers have the address 114 and are labelled as slave controls.

Each controller has to have an individual address. No address may be repeated.

Master control has only got the address 113.

Slave control begins with adress 114 and ends with adress120.


The actual control address is shown on the display of the controlbox SM7.

S	y	s	t	e	m		r	e	a	d	y					
A	d	r	:	1	1	3	-	M	a	s	t	e	r			

The control address can be changed by using the programme „LN-SM7-Configuration Tool“.

Cross-link a Slave control with a Master control

- Switch off both controls.
- Plug one end of the LAN cable in the Output bush of the Master control and the other end in the Input bush of the Slave control.
- Switch on both controls again.

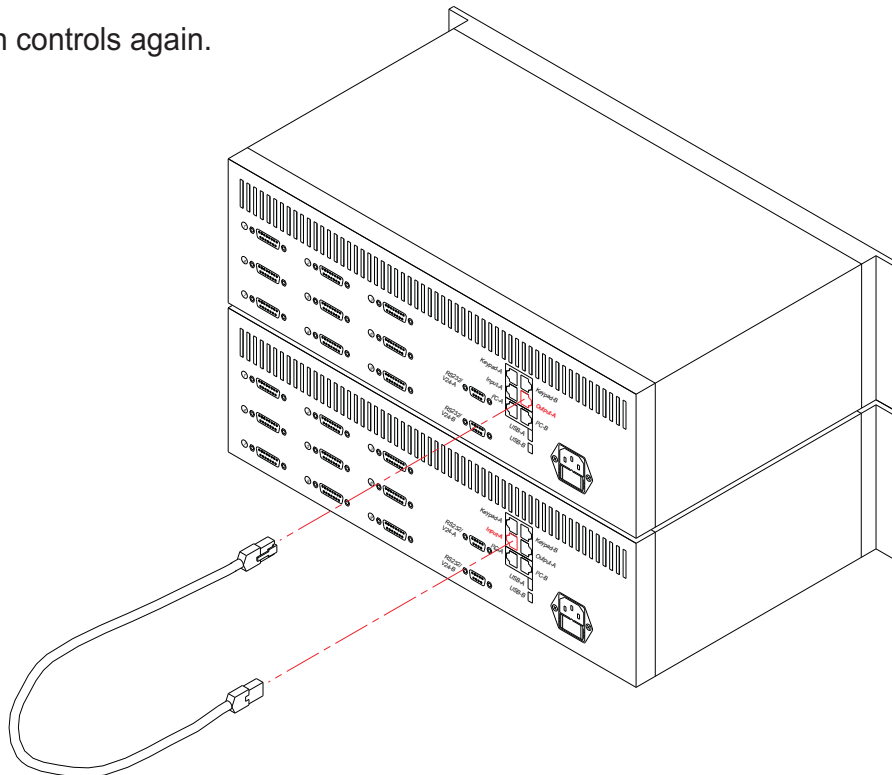

Master = The first connected controls in the network-chain, set to address 113 and attached to the keypad SM7.

Slave = All following controls in the network-chain. The slave control starts at address 114. One master control can be networked with up to seven slave controls. No slave control can be connected with a keypad.

LN-SM7 ConfigurationTool

Each control SM7 has fixed physical addresses according to the number of motor drivers. These physical addresses are related to virtual addresses. The configuration of the devices follows automatically. Altogether eight controls can be cross-linked. This complies with maximum 72 motor drivers / motor-driven axes.

The programme “*LN-SM7 ConfigurationsTool*” is used for the configuration of the virtual addresses. A condition for using installed at Microsoft .NET Framework 4 Extended or higher. Check under System control (Systemsteuerung)Programme (Programme).

Should a further control be connected with a control SM7, the address on the additional control must be changed. The first control which is connected with the keypad has got the label Master control and the address 113. The subsequent controls have got the label Slave control and the addresses from 114 to 120.

No address may be repeated.


The control address can be changed by using the “LN-SM7 ConfigurationTool”

Changing address

- Connect the controller to the computer over USB and switch on the control SM7




When addressing the controller to the computer, the control may not be connected with another control.

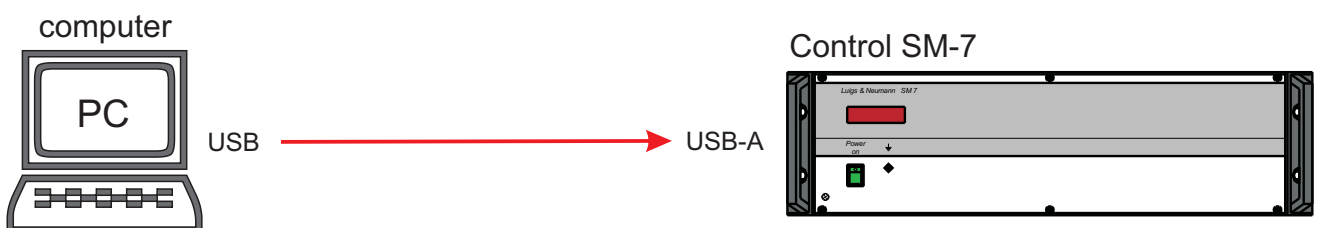


Setting the control address

Start “LN-SM7-ConfigurationToll“

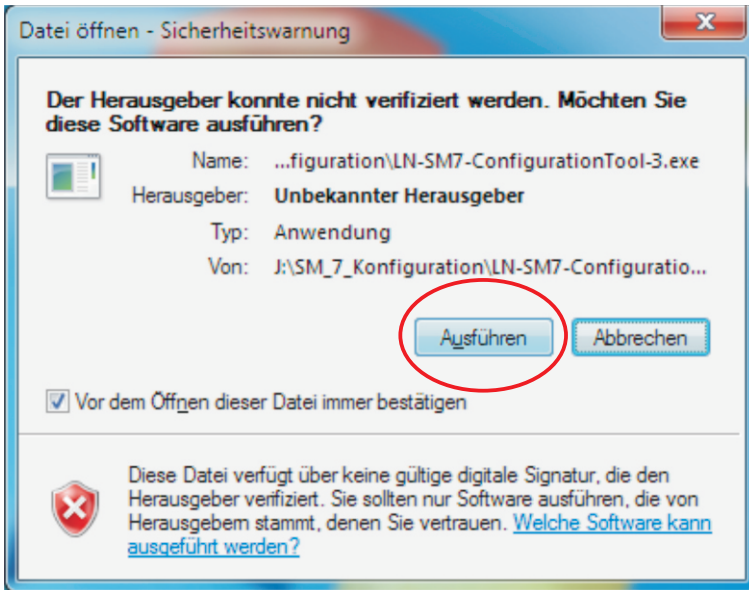

Click on “Ausführen“

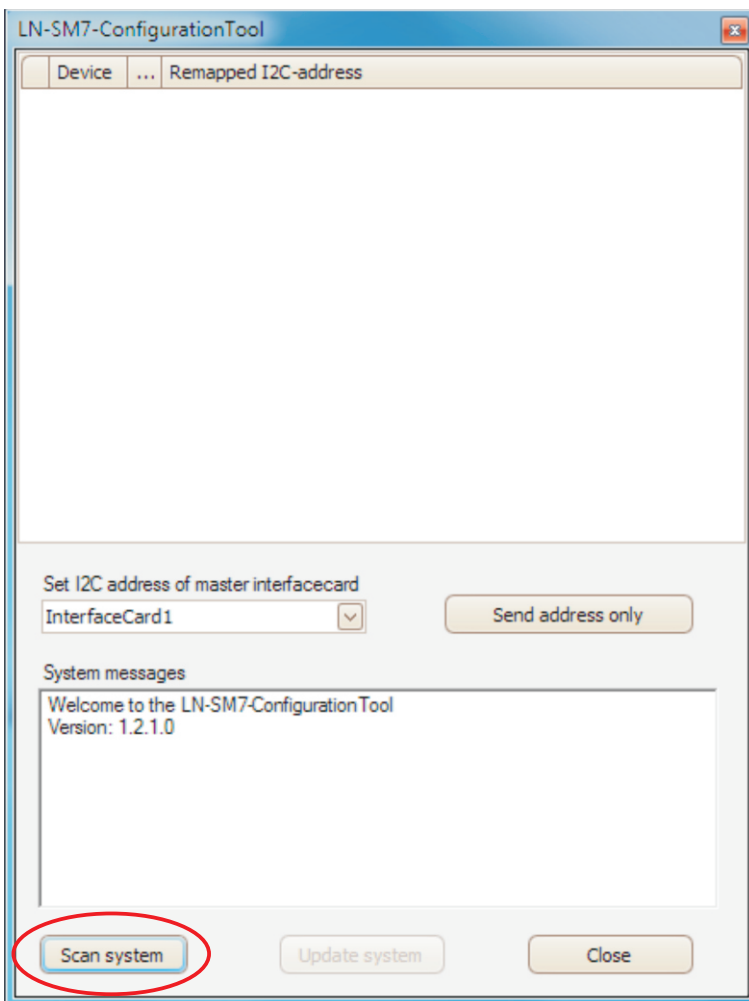


Click on “Scan system“

Setting the control address

The result is shown:

Example: Control SM7 with 9 motor-driver and the address 113

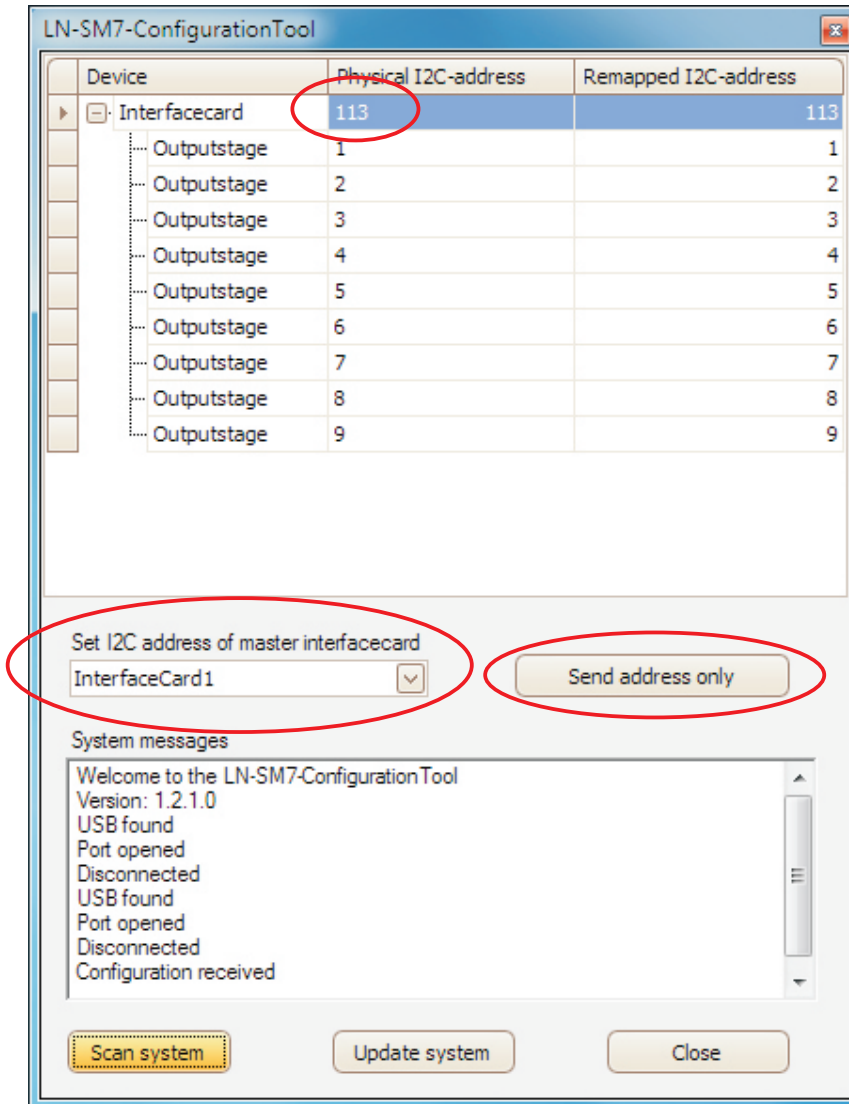

Relation of addresses:

InterfaceCard 1 = **113** (Master)
InterfaceCard 2 = **114** (Slave)
InterfaceCard 3 = **115** (Slave)
InterfaceCard 4 = **116** (Slave)
InterfaceCard 5 = **117** (Slave)
InterfaceCard 6 = **118** (Slave)
InterfaceCard 7 = **119** (Slave)
InterfaceCard 8 = **120** (Slave)

The actual address is shown in the line "Interfacecard".

In the line "**Set I2C address of master interfacecard**" the desired address is set and confirmed by clicking "**Send address only**".


After the restart of the control SM7, the new address is taken over.

Example: Address the control on 114 and configuration the motor connections

Example:

Two controllers with eight motor-drivers should be installed for a work station with four 4-axes manipulator units.

The first control which is connected with the keypad Sm-7 must have the address 113 and the second control the address 114.


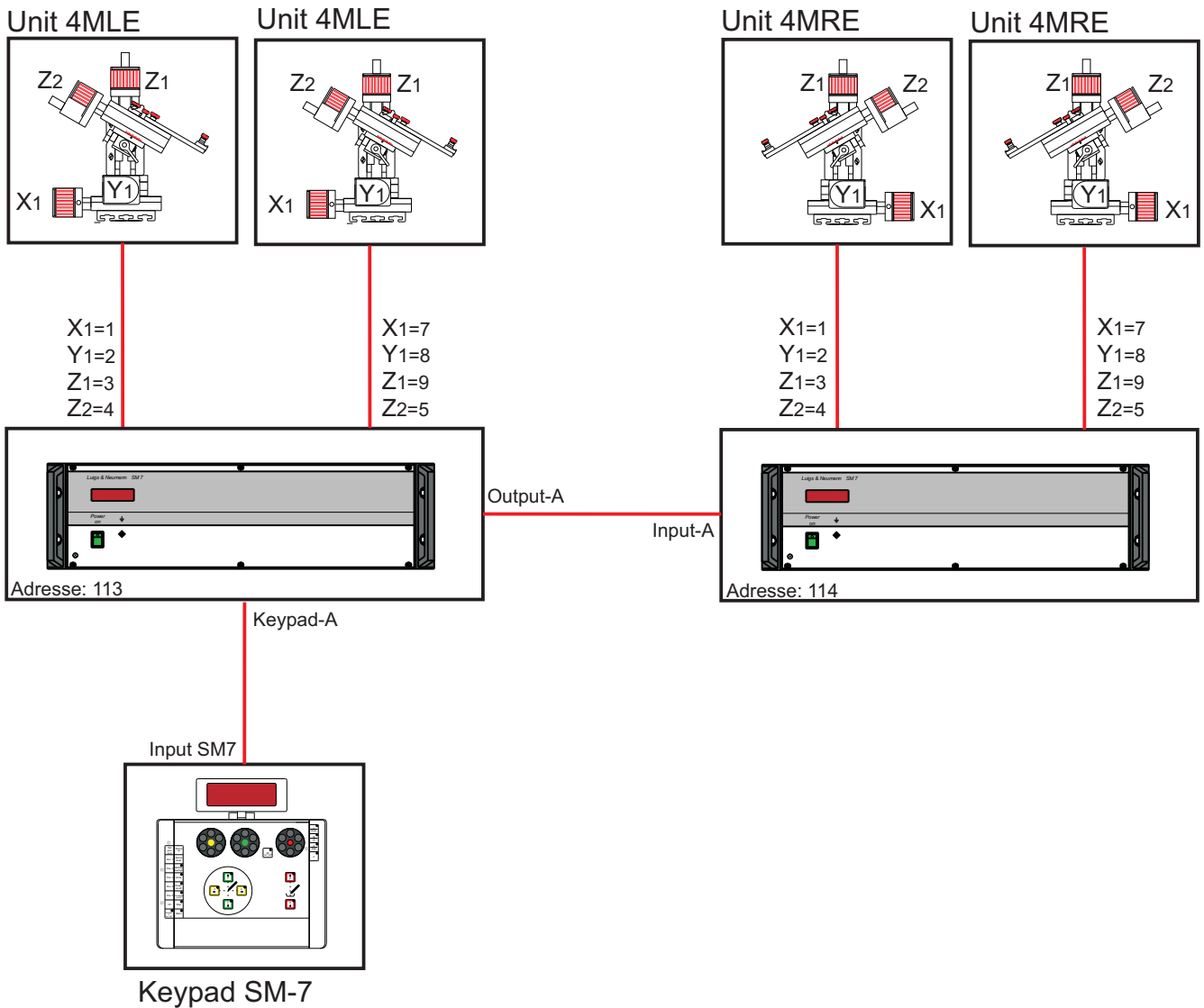


Example: Address the Slave-control on 114

Connect the controller to the computer over USB.



Select and open the programme “LN-SM7-ConfigurationTool”

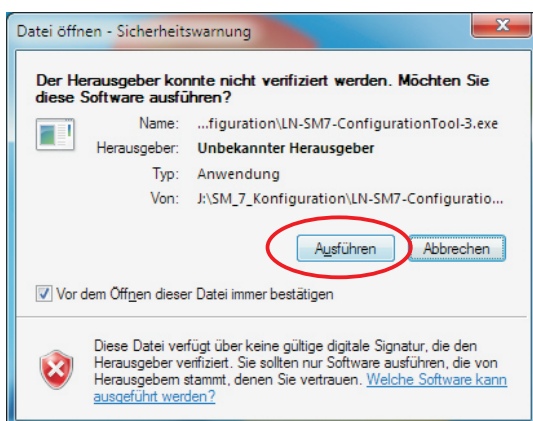
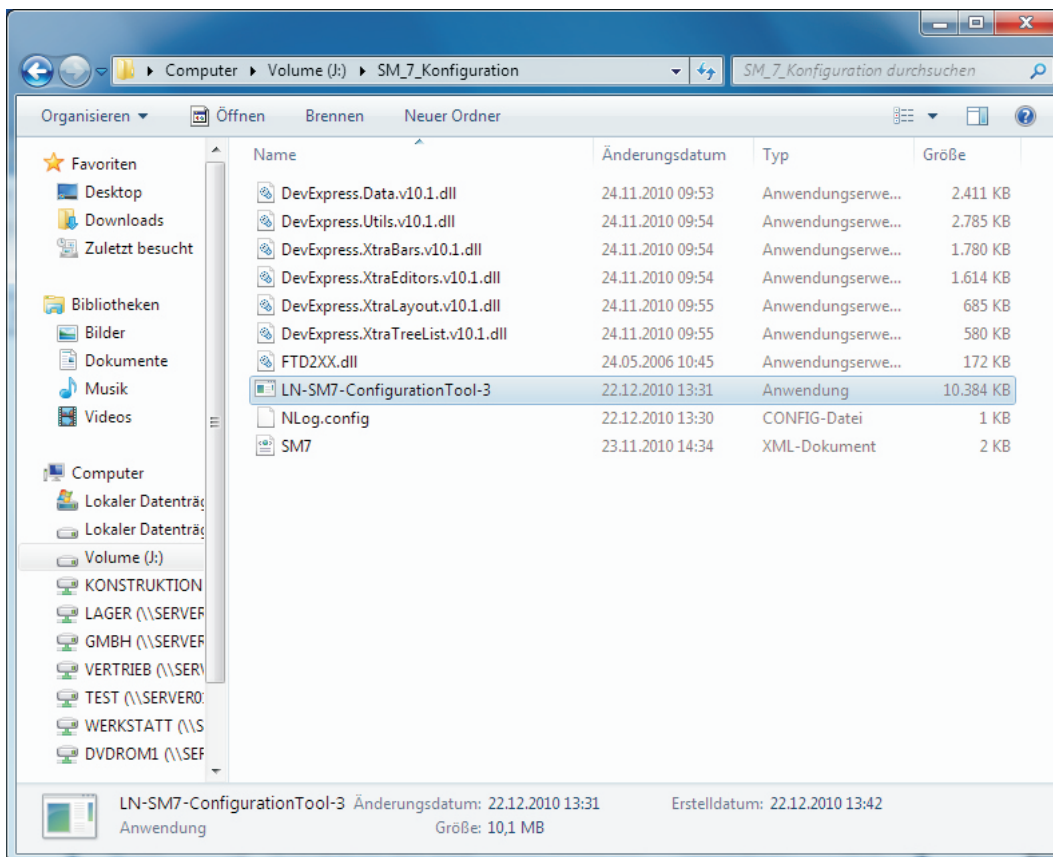


Click on “Ausführen”



Example: Address the Slave-control on 114

Set the address 114 for the second control:

Additionally, in the line „ **Set I2C address of master interfacecard**“ select InterfaceCard2

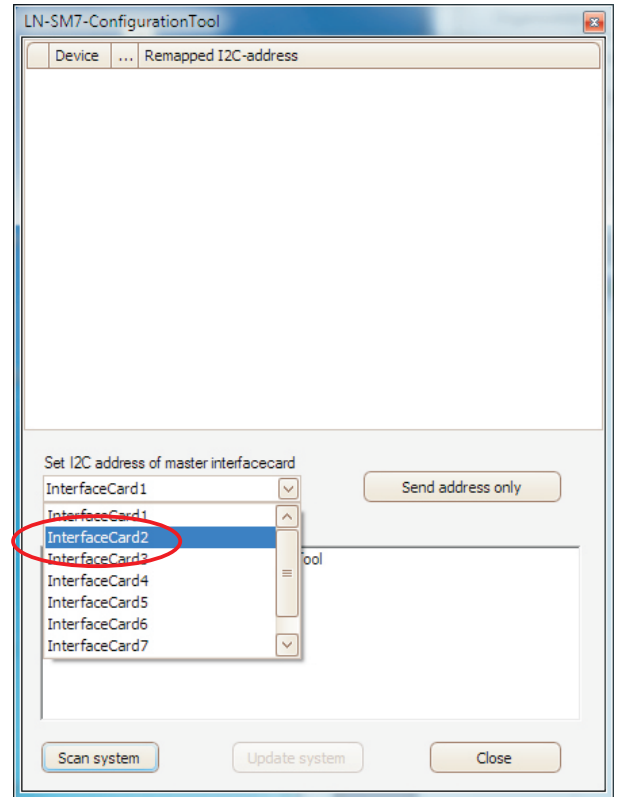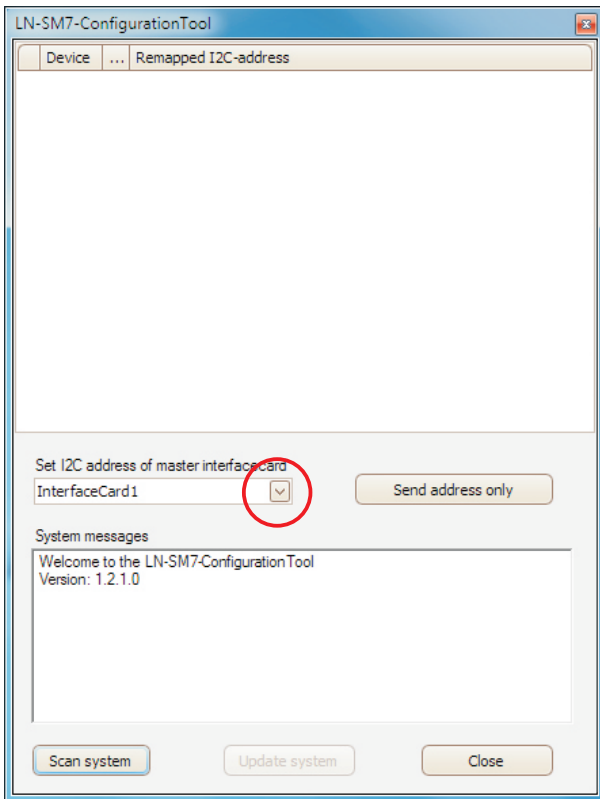


and click on **“Send address only”**.

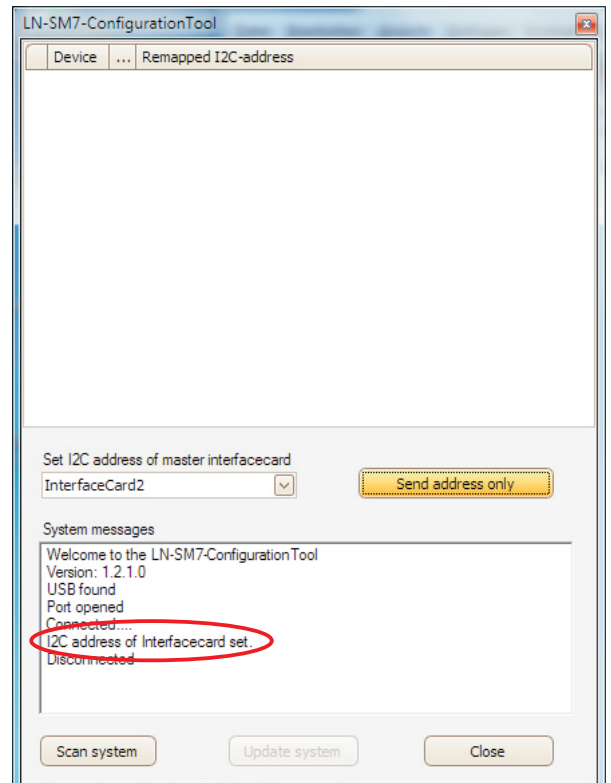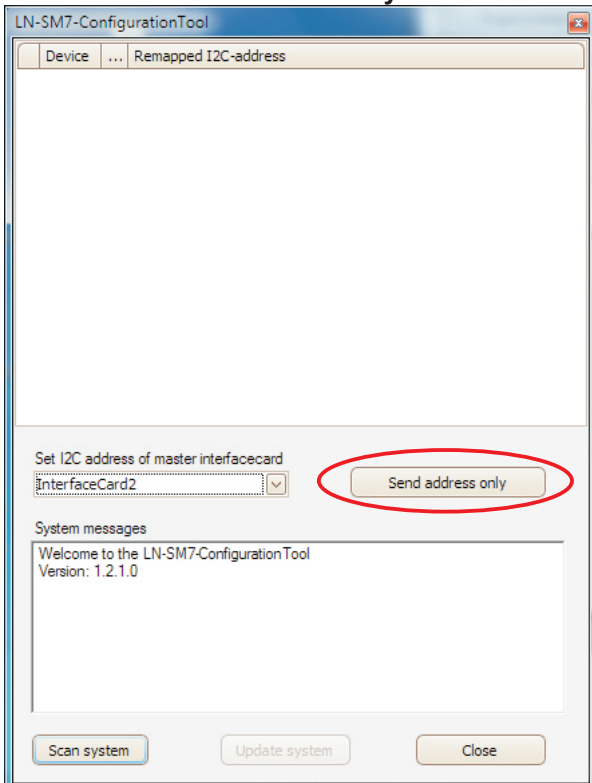


In the line **“System messages”** the confirmation of changes is shown: ***I2C address of Interfacecard set.***

Example: Configuration motor connections

- Close the “LN-SM7-ConfigurationTool” and switch off both controls.
- Plug one end of the LAN cable in the **Output-A** socket of the master control and the other end in the **Input-A** socket of the slave control.

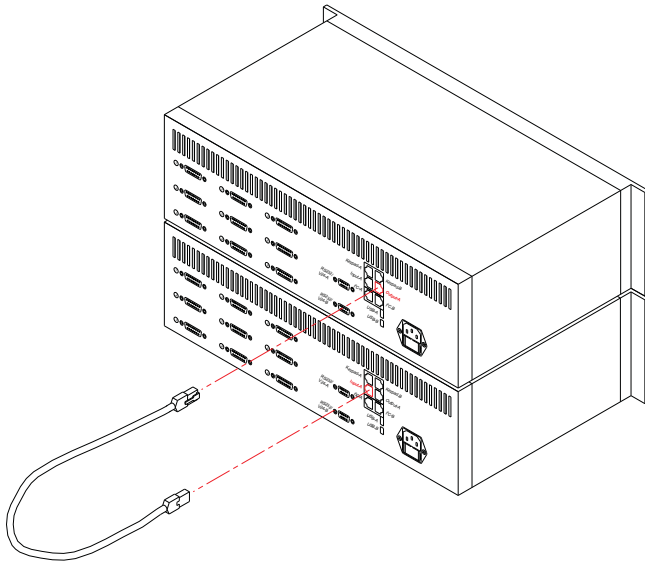



Connect the controller to the computer over USB

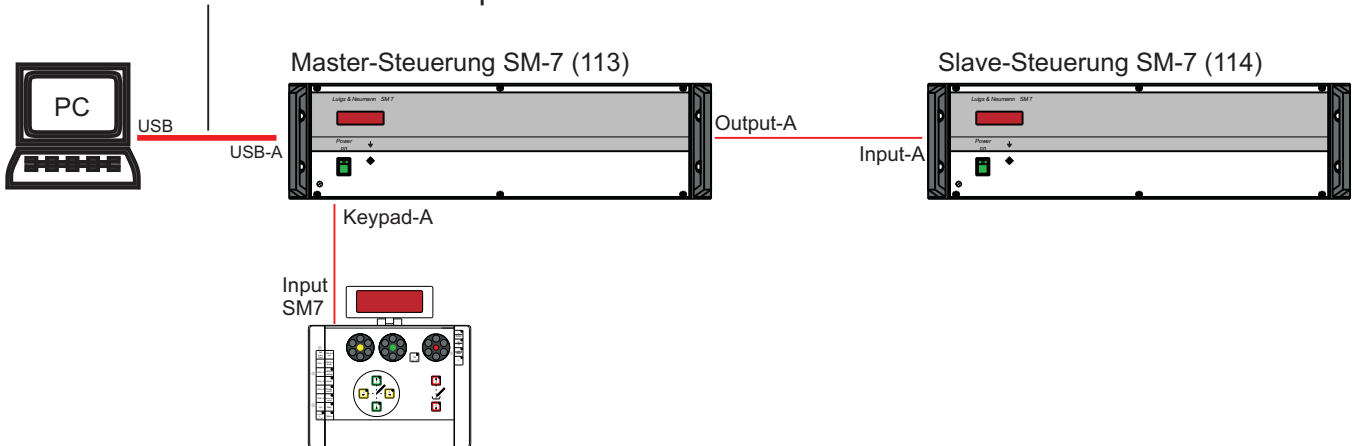

Switch on both controls and start “LN-SM7-ConfigurationTool”.

Example: Configuration motor connections

Click **“Scan system”**

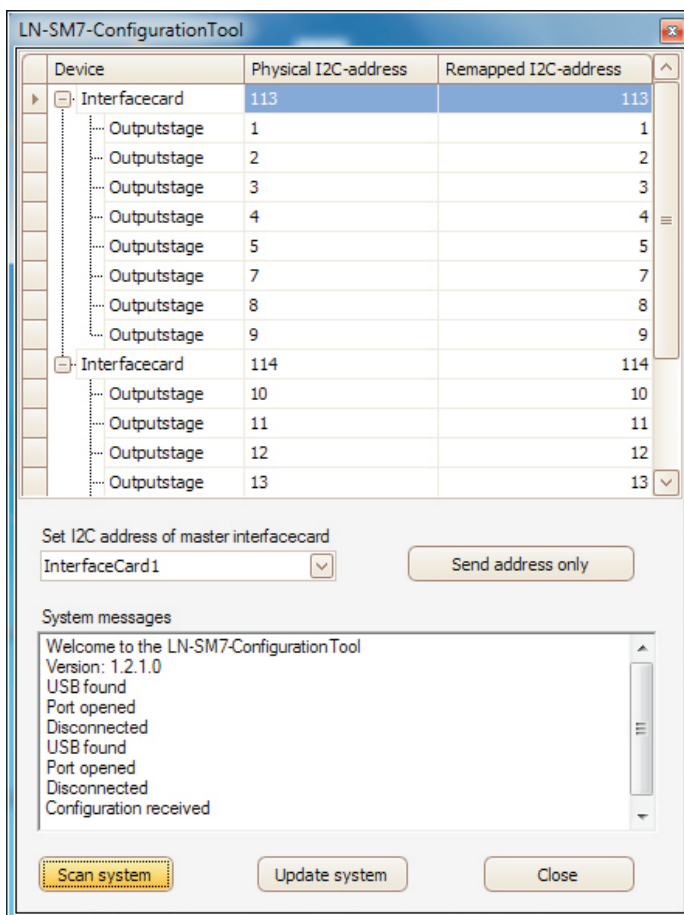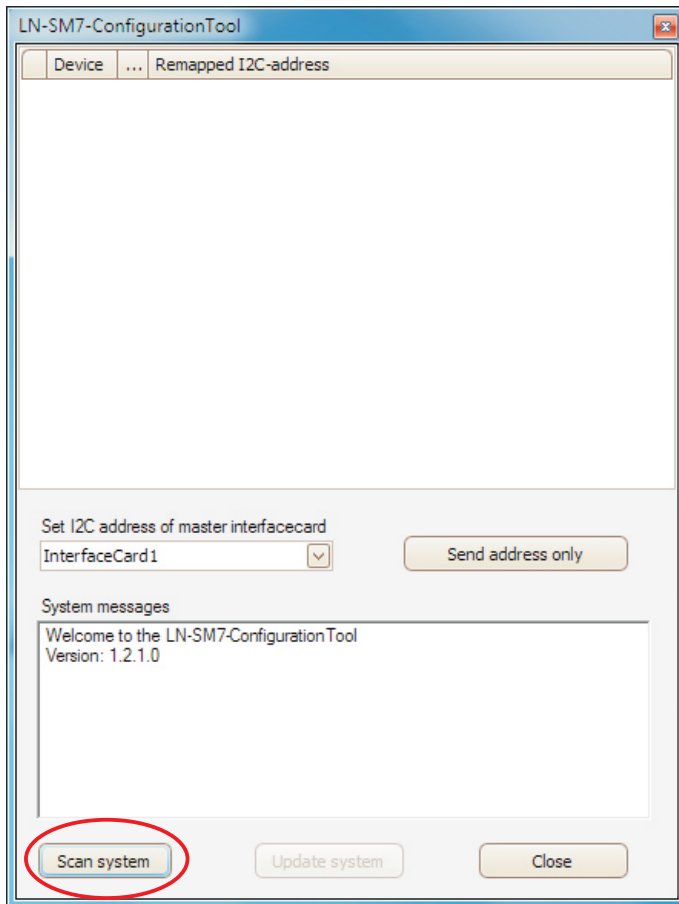

The configuration of the motor connections is shown.


Example: Configuration motor connections

The keypad SM-7 should show the following virtual configuration.
 For that the single motor connections in the control SM7 must be set up with the programme "LN-SM7-ConfigurationTool".

Keypad SM7: Display position-No.

Dev 1	X 1	Y 2	Z 3
Dev 2	X 4	Y 5	Z 6
Dev 3	X 7	Y 8	Z 9
Dev 4	X 10	Y 11	Z 12
Dev 5	X 13	Y 14	Z 15
Dev 6	X 16	Y 17	Z 18
Dev 7	X 19	Y 20	Z 21
Dev 8	X 22	Y 23	Z 24


Example:

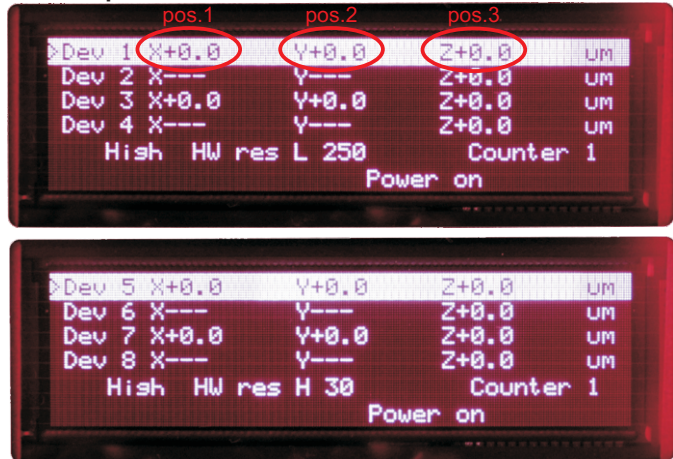


Virtual co-ordination of display position

	Axes	Remapped I2C-adress Display-position Keypad SM-7	Physical I2C-adress Motor connection Control SM-7
1. Unit MLE	X1	1	1
	Y1	2	2
	Z1	3	3
	Z2	6	4
2. Unit MLE	X1	7	7
	Y1	8	8
	Z1	9	9
	Z2	12	5
3. Unir MRE	X1	13	10
	Y1	14	11
	Z1	15	12
	Z2	18	13
4. Unit MRE	X1	19	16
	Y1	20	17
	Z1	21	18
	Z2	24	14

Physical co-ordination of motor connection
 1. control: Master

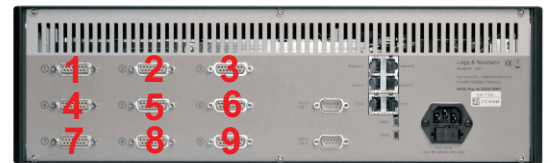


Physical co-ordination of motor connection
 2. control: Slave

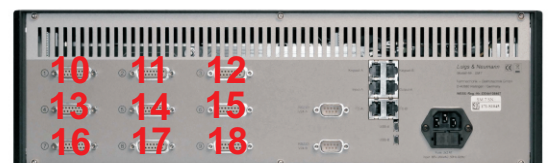

Example: Configuration motor connections

Enter the desired virtual co-ordination of the single axes in the column “Remapped I2C-address”

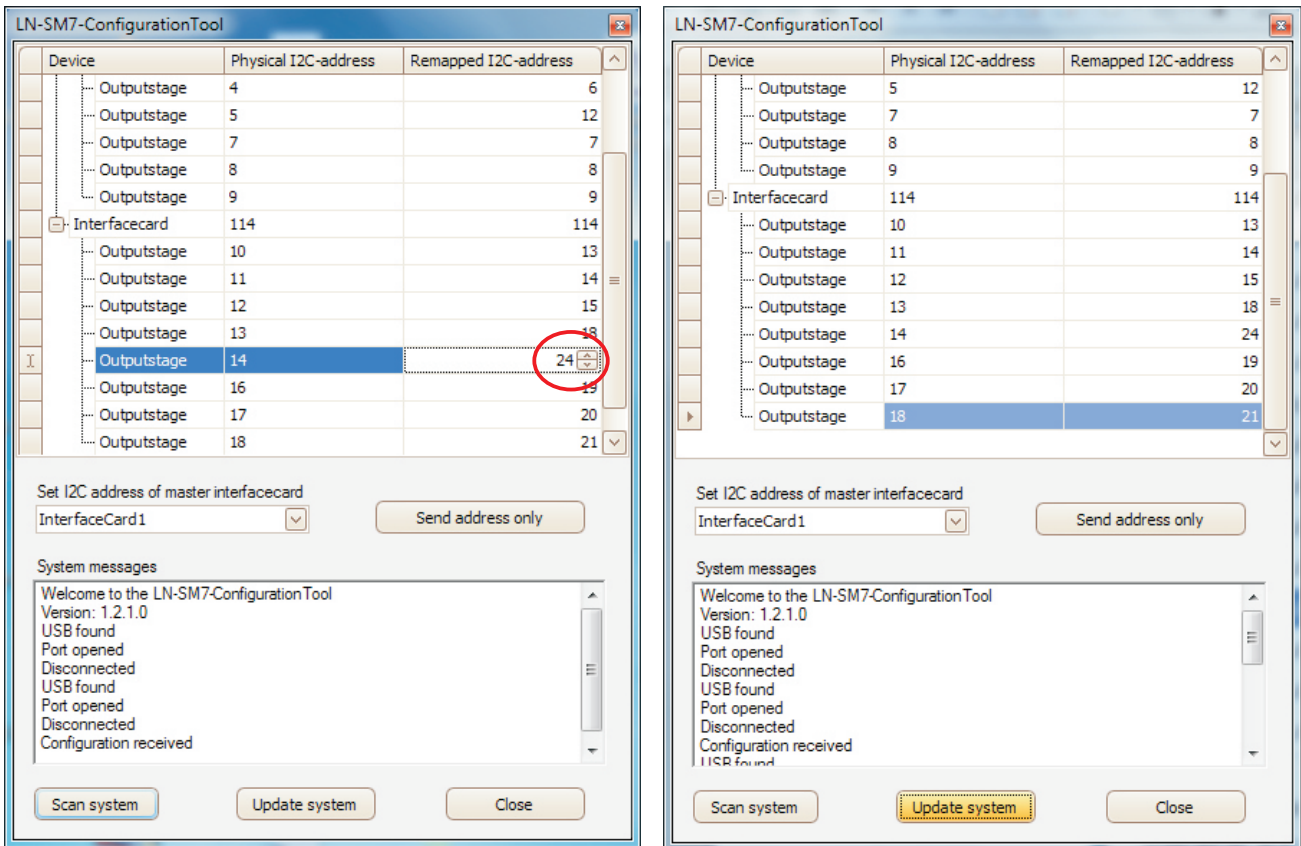


Click “Update system”

Switch both controls off and then on again.

The desired axes configuration is shown on the display of the keypad.



Appendix - B

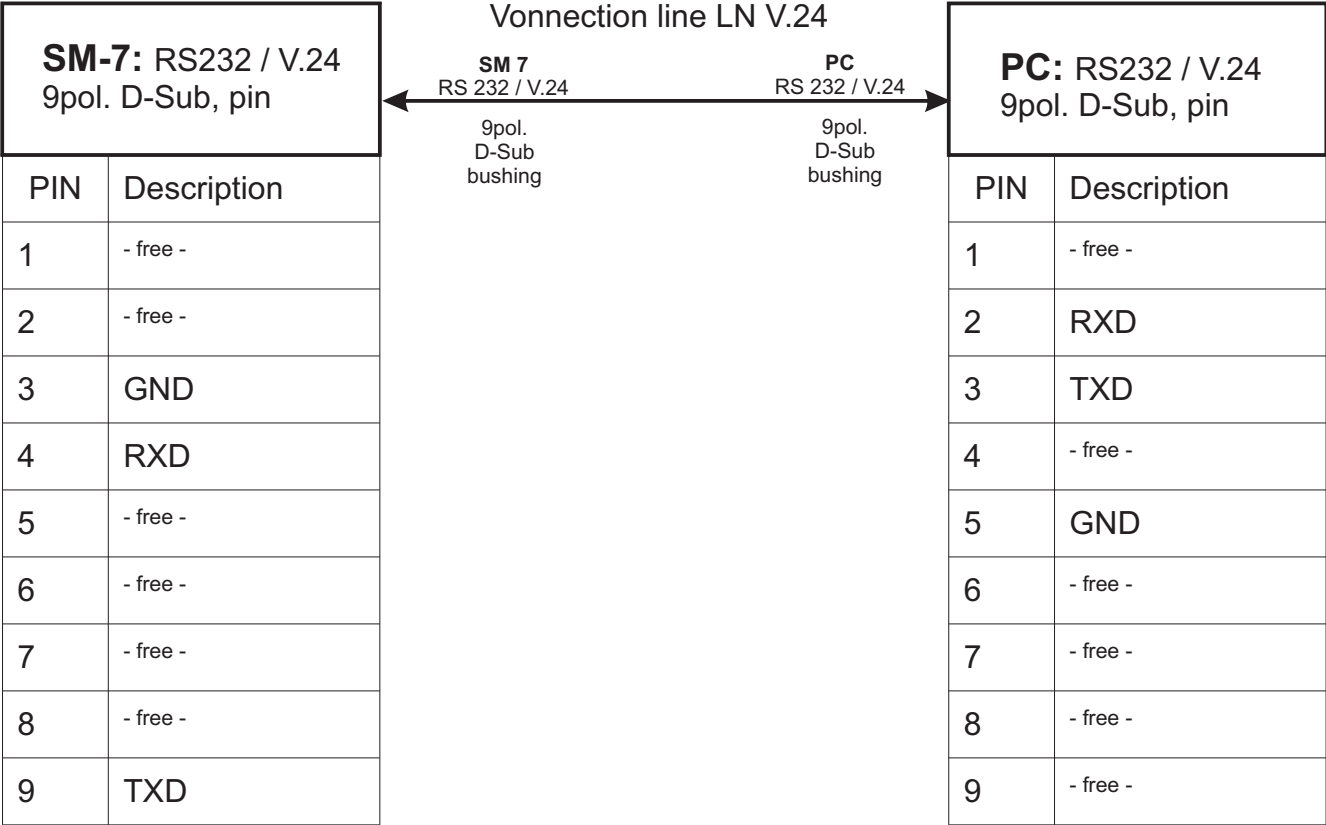

Technical Data

Casing	: 19" / 3HE, 255,5mm deep
Exterior dimensions	: 449mm x 133mm x 255,5mm
Weight	: 6,5kg
Input voltage	: 85V – 264V AC, (47-63Hz)
Fuse	: 2x T2A
Performance	: 150W
Current	: max. 1,3A bei 100VAC / 0,65A bei 200VAC
Inrush current (100V/200V AC)	: 14 / 28A
Ambient temperature	: 0°C – 40°C
Relative humidity	: 90%

Applicable law:
EMV Richtlinien 04/108/EG
Niederspannungsrichtlinie 72/23/EGW
Ergänzt durch 93/68/EGW